Experimental Procedure Preparation of Methyl N-tert-Butoxycarbonyl L-Alanyl-L-Phenylalaninate

- 1. Place 0.63 g of methyl L-phenylalaninate hydrochloride in a 150-mL Erlenmeyer flask.
- 2. Add 10 mL of dimethylformamide and 0.3 mL of N-methylmorpholine to the flask, swirl the flask to mix the contents, and place the flask in an ice-water bath. Label the flask as Solution A.
- 3. Add 0.50 g of N-tert-butoxycarbonyl-L-alanine in a 100-mL round-bottom flask containing a one inche stirbar. Add 10 mL of dimethylformamide, and then using a 1-mL syringe transfer 0.3 mL of N-methylmorpholine to the flask. Place the flask in an icewater bath and stir the solution for 5 min.
- 4. Using a 1-mL syringe add 0.4 mL of isobutyl chloroformate dropwise to the cooled solution in the round-bottom flask and continue stirring the mixture in the ice-water bath for 5–10 min.
- 5. Transfer Solution A into the round-bottom flask and continue to stir the reaction with cooling in the ice-water bath for 45 min.
- 6. Add 20 mL of water to the reaction mixture and transfer it to a separatory funnel. Rinse the round-bottom flask with a 30-mL portion of diethyl ether and transfer the rinse to the separatory funnel.
- 7. Shake the funnel gently, venting as necessary. Separate the layers and wash the organic layer with two 25-mL portions of 1 M HCl, a 25-mL portion of saturated sodium bicarbonate, and a 25-mL portion of brine.
- 8. Transfer the organic layer to a 125-mL Erlenmeyer flask and add several spatula-tips full of anhydrous sodium sulfate.
- 9. Decant the solution into a 150 mL beaker and air dry the sample until the ether is almost gone. Add 2-3 mL ether back into the beaker and swirl until the solid has dissolved.
- 10. Add approximately 15 mL of hexanes to the flask and place the flask in an ice-water bath for about 15 minutes. Use a spatula to scrape all the solid from the bottom of the flask.
- 11. Isolate the crystals by vacuum filtration and air-dry them.
- 12. Weigh the recrystallized product, determine its melting point, and obtain an IR spectrum.